

# **STANDARD OPERATING PROCEDURES**

5. **Check for sufficient printer paper**, if red ink is showing, reload with new roll.
6. **Close chamber door** using touch screen display by pushing "**CLOSE DOOR**". Maintain pressure on display until the door is closed.
7. Choose main menu, and then **select appropriate cycle** option(s).
8. Daily air removal tests should be performed each day prior to running a load. If the autoclave has not been used in several days, press "**CYCLE SELECT**" select "**TEST CYCLES**" and select "**DART WARMUP**" cycle to bring chamber up to operating temperature in preparation for the DART/Bowie Dick test cycle.
9. After "**DART WARMUP**

SOP# 1150.1  
Amsco Model #85 Century Bulk Sterilizers  
Effective 7/13